




User manual

- 2 -

- 1 -

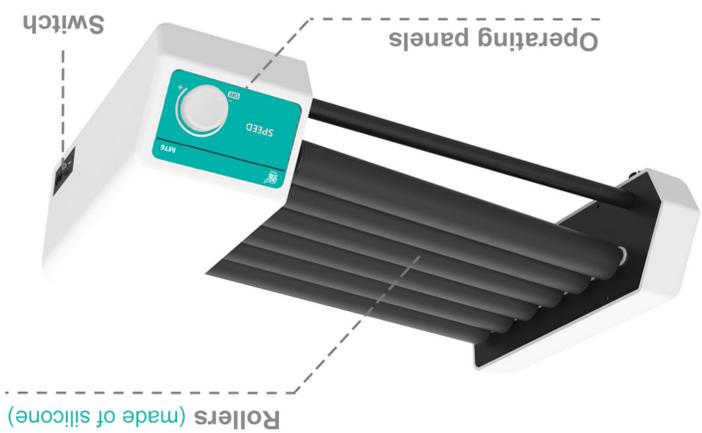

II: Product overview

This mixers are designed and manufactured according to international advanced design concepts and technologies. The mixers are highly efficient and gentle and are suitable for different experiments such as blood analysis and sample mixing.

III: Product introduction

I: Product brief

IV: Product data

Model	MT6
Products Name	Roller mixer (High-low axis)
Image	
Type	Rotating shaking
Rollers	6 rollers
Max Capacity	12*50ml Centrifuge tube / 0.6kg
Control mode	Digital Key
Speed	20-80 rpm
Timing	/
SPECS	φ6-34mm
Power	25W
Voltage	AC100-220V
Dimension(mm)	440*255*105

V: Correct usage

/// Basic usage

Application

> Suitable for mixing operations in different applications: electron microscopy analysis, blood cell suspensions, homogeneous solid blood and plasma, tissue specimens, washing precipitates, disintegration sheet analysis, extraction process, dissolution slow reaction, rotary sealing, etc.

Usage

> Widely used in scientific research institutions, medical schools, health care and other units, molecular biology, virology, microbiology, pathology, immunology and other laboratories.

- 3 -

Delivery list


> Host > Power cable > User manual



+



+

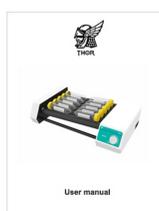

- 5 -

Steps for usage

- > Plug in the power supply and turn on the switch;
- > Place the sample tubes on the roller;
- > Turn the knob to adjust the rotational speed;
- > Turn the knob to the left to turn off the device.

▲ Instrument using environment

- > Indoor use
- > Altitude: ≤2000m
- > The instrument is suitable for ambient temperature range of +5°C~+40°C
- > Instrument operation suitable for relative humidity range of ≤80%
- > No vibrations and airflow affecting performance are present
- > There are no conductive dust, explosive gases and corrosive gases in the surrounding air
- > Keep a safety distance of 10cm between the front and back left and right when the instrument is working

▲ Safety Precautions

- > It is strictly forbidden to unplug the power connector and toggle the switching power button when your hands are stained with liquid!
- > It is strictly forbidden to unplug the power plug when the instrument is live!
- > It is strictly forbidden to maintain and scrub the instrument when it is powered on!
- > It is strictly forbidden to install the instrument on an uneven, shaky and shaky work surface!

- 4 -

Product conformity certificate

Product Name : Roller mixer

Model number : MT6

Production date : _____

- 6 -